VMS-D
VMS-A

VMS-A

VMS-D

Betriebsanleitung	DE	5
Operating instructions	EN	14
Mode d'emploi	FR	23
Istruzioni per l'uso	IT	33

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Warranty EN

In accordance with **VWR** Terms and Conditions of Sale, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the device direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance

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Safety instructions

Do not use the device in explosive atmospheres, it is not EX-protected.

With substances capable of forming an explosive mixture, appropriate safety measures must be applied, e.g. working under a fume hood.

To avoid body injury and property damage, observe the relevant safety and accident prevention measures when processing hazardous materials.

- Set up the device in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.

Attention – Magnetism! Effects of the magnetic field have to be taken into account (e.g. data storage media, cardiac pacemakers ...).

- Ensure that the power cord set / temperature sensor cable does not touch the heating plate.
- **Version VMS- D:** Examine regularly (we recommend every 4 weeks) the function of the safety temperature limiter. For this the safety temperature must be changed around at least 1/4 revolution of the value already stopped and be accomplished the attitude as under “adjusting the safety temperature delimitation SAFE temp” described. By the exceeding of the adjusted safety temperature (e.g. rotary button “temp” posed on right end stop) you can examine the function. The “temp” pilot lamp must expire and the heating must be lasting switched off.
- When using PTFE-coated magnetic bars, the following has to be noted: Chemical reactions of PTFE occur in contact with molten or solute alkali metals and alkaline earth metals, as well as with fine powders of metals in groups 2 and 3 of the

periodic system at temperatures above 300 °C – 400 °C. Only elementary fluorine, chlorotrifluoride and alkali metals attack it; halogenated hydrocarbons have a reversible swelling effect.

(Source: Römpps Chemie-Lexikon and “Ullmann”, Volume 19)

Risk of burns! Exercise caution when touching parts of the housing and the heating plate. The heating plate can reach dangerous temperatures. Pay attention to the residual heat on the heating plate after switching off the stirrer.

The device may only be transported when the heating plate has cooled down.

- Please consider any possible contaminations and unwanted chemical reactions.
- Before placing the device in service, please ensure that the turn dial for adjusting the speed is against the left stop, since the device will begin to run at the last speed that was set.
- The device will automatically restart following any interruption to the power supply.
- The device may not be transported by using an H 16 V support rod or the power cable.
- Clicking noises may occur as a result of material expansion during heating and cooling.
- **Warning!** The safety temperature must be set in accordance with EN 61010-2-010 Chapter “Requirements for devices containing or using flammable liquids”.
 - The surface temperature of the flammable medium that is exposed to air may not exceed its flash point. A danger usually arises if a medium is heated in open vessels.
 - The surface temperature of the heating device (e.g. the mounting plate) may not exceed the value of $(t - 25)$ °C (= set value of the safety circuit) on the surface of the flammable medium and in contact with air, whereby t is the fire point of the liquid. A danger usually arises if a medium is heated in glass vessels (glass breakage).

If a setting made by the user (medium temperature or safety temperature) could bring a flammable medium into a state in which the conditions mentioned above could be exceeded, additional measures must be introduced that will protect the user from this danger.

- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- The device may only be opened by qualified and VWR approved experts.
- Safe operation is guaranteed only with the use of original VWR accessories.

Intended use

Use

- The VWR VMS- A or VMS- D is a stirring instrument with heating function. The instrument is suitable for tempering substances which are placed on the heating plate in vessels. By means of a built-in stirring drive and with the help of a magnetic bar within the vessel, the substances can be stirred at the same time. The mixing intensity is dependent on the motor speed and on the size of the magnetic bar. The VMS- D also has a contact thermometer connection and an adjustable safety circuit.

Area of use

- Indoor environments similar to that a laboratory of research, teaching, trade or industry area.
- The safety of the user cannot be guaranteed:
 - if the device is operated with accessories that are not supplied or recommended by the manufacturer,
 - if the device is operated improperly or contrary to the manufacture’s specifications,
 - if the device or the printed circuit board are modified by third parties.

Contents

	Page
EU Declaration of conformity	3
Warranty	4
Safety instructions	14
Intended use	15
Unpacking	16
Commissioning	16
Starting	16
Function stirring	17
Function heating	17
Limiter of safety temperature	18
Adjusting the safety temperature	
limit “Safe Temp” (Version - D)	18
Controlling the medium temperature limit using a contact thermometer	19
Error messages	20
Mounting support rod H 16 V	21
Maintenance and cleaning	21
Accessories	21
Technical Data	22

Unpacking

Unpacking

- Please unpack the device carefully
- In the case of any damage a detailed report must be sent immediately (post, rail or forwarder)

Scope of delivery

- VMS- A or VMS- D
- Operating instructions

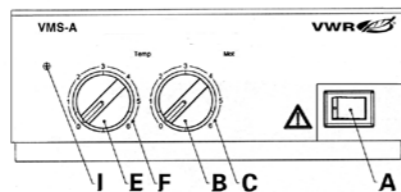
Commissioning

Check whether the voltage specified on the type plate matches the mains voltage available. The power socket used must be earthed (protective earth conductor contact). Before turning on the machine, adjust the turn dial for controlling the speed so that it is against the left stop. If these conditions are met, the device is ready to operate after plugging in the mains plug. If these procedures are not followed, safe operation cannot be guaranteed and/or the equipment may be damaged. Observe the ambient conditions (temperature, humidity, etc.) listed under "Technical Data".


Starting

The heating stirring features can be turned on and off by pressing the unit switch **A**.

Version VMS- A: The ready for operation and turned on states of the device are indicated by the green lamp in the unit switch lighting up.

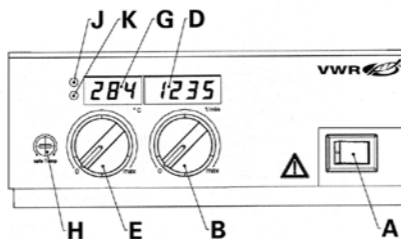Version VMS- D: When you switch on the stirrer it will run a selftest and all the LED segments of the displays will light up for a few seconds. After this, the LED's (**D**; **G**) light up to indicate that the unit is powered up and ready for operation.

Function stirring

The stirring rod in the glass beaker with a maximum length of 60 mm is driven by a permanent magnet located directly under the placing and heating surface. It allows magnetic field lines to pass through it. The permanent magnet is fastened directly onto the drive shaft of the motor.

The actual speed depends on the load and voltage. Note that fluctuations in the in the mains power voltage within the allowable tolerance and changes in viscosity of the materials to be stirred resulting from processes they undergo may also cause minor fluctuations in the speed.

Version VMS- A: The speed of the motor can be adjusted on the "Mot" dial (**B**) with a scale from 0 to 6 (**C**). The scale values from 0 to 6 on this dial correspond roughly to speeds ranging from 0 to 2000 rpm. The maximum speed of 2000 rpm is reached when the dial switch is turned all the way to the right and there is no load on the motor.

Version VMS- D: Set the motor speed using the "Mot" knob (**B**); the knob has a graduated scale from to maximum rpm.

The motor speed is shown on the LED display (**D**). When the motor is stopped the speed "0" will be displayed. When the motor is operating, the display will show the current speed in the range 100 to 2000 rpm.

Function heating

The temperature of the heating plate of the instrument is kept constant by a control circuit. In addition, it is monitored by a safety circuit. The required temperature sensor (a thermo-couple) is installed in the heating plate. The **VMS- A** unit has a stainless steel surface plate with a heating output of 400 Watt.

VMS- D units have a enamel-coated stainless steel surface plate with a heating output of 500 Watt.

Version VMS- A: The temperature of the heating plate can be adjusted on the „Temp“ turn dial (**E**) with a corresponding scale from 0 to 6 (**F**). The scale values from 0 to 6 correspond approximately to area temperatures up to a max. of 320 °C. When the setting is at 0, the device does not heat the medium up at normal room temperature.

During the heating phase, the signal light (**I**) lights up orange (orange = power is being supplied to the heating plate; signal light off = no power being supplied)

Version VMS- D: The temperature of the heating plate can be adjusted on the „Temp“ turn dial. When you change the knob setting, the LED display (G) will show the SETPOINT temperature and the warning light (K) will light up. After approx. 3 seconds the LED display will switch back to showing the ACTUAL temperature and the warning light (K) will switch off.

The temperature indication now constantly changes between actual temperature (25 sec) and target temperature (5 sec).

When you select the SETPOINT temperature and as the unit heats up, the warning light (K) will remain off and the LED display (G) will show the ACTUAL temperature of the surface plate in the range from 0 °C to + 320 °C.

The red warning light (J) will light up while heating is in progress.

(Red warning light ON = surface plate heating power supply is ON; red warning light OFF = surface plate heating power is OFF).

In position "0" the device does not heat at room temperature. If, however, the temperature of the heating plate is higher than 50 °C, the LED display (G) flashes to indicate "Hot".

PLEASE NOTE !

The heating plate of the version **VMS- D** only heats up, if the contact plug has been plugged (the contact plug is undetachably fitted to the instrument), or if VT 4 or an contact thermometer has been connected.

Limiter of safety temperature

The safety temperature limiter makes it possible to adjust the maximum temperature of the heating plate (safety temperature).

This ensures that the temperature of heating plate does not exceed the flame point of the material. The value to be set must be determined by trial and error and should generally fall about 20 K below the flame point of the medium.

VMS- A devices have a safety circuit with a fixed setting that limits the temperature of a heating plate to 400 °C.

VMS- D devices have an adjustable safety circuit. The maximum temperature of the heating plate can be adjusted within a range from 100 °C to max. 400 °C.

Adjusting the safety temperature limit "Safe Temp" (Version - D)

Version VMS- D: The "Safe Temp" set screw (H) is to the left of the "Temp" knob (E) on the front operator panel.

You can use the „Safe Temp“ set screw (H) of the safety temperature limiter to adjust the desired safety temperature. If the desired safety

temperature is exceeded by a higher adjusted target temperature, or if there is a defect in the instrument controller, the heating turns off permanently in this state and the stirring function continues to run at the speed that was set prior to the malfunction.

- Using a straight-edge screwdriver, turn the "Safe Temp" adjustment screw (H) as far as it will go clockwise (to the right).
- Do not turn the screw beyond the natural right or left limit of rotation, as this will destroy the potentiometer.
- Use the "Temp" turn dial (E) to adjust the set-point temperature to desired safety temperature ("Safe Temp").
- Turn the "Safe Temp" set screw (H) slowly to the left until the heating function switches off and the signal light (I or J) goes off. Wait for about one more minute to ensure that the signal light (F) going off does not merely indicate the end of a switching cycle of the heater. If the signal light (I or J) turns on again within this time, you must repeat the procedure, slowly turning the "Safe Temp" (H) adjustment screw to the left again. When the signal light (I or J) does not light up again within one minute, the safety temperature is correctly set.
- After this, turn set "Safe Temp" adjustment screw (H) back slightly to the right (clockwise).
- After you have turned the unit switch (A) off and back on again, the device is ready for operation.

NOTE: When the "Safe Temp" (H) adjustment screw is set to the minimum level (as far as it will go to the left) the heating function is turned off.

Triggering of the safety circuit is shown in the following way:

Version VMS- A: The warning light (I) will flash two times per second. (Note: regular flashing indicates that the safety circuit has been triggered; this warning light remains permanently lit to indicate that a standard heating cycle is in progress).

Version VMS- D: The letter "-5-" will be shown on the LED display (G).

Controlling the medium temperature limit using a contact thermometer

The preferable method for controlling the average temperature is with contact thermometer. After the set point temperature has been adjusted, this results in a short heating-up time, practically no temperature drift and only minor fluctuation in temperature.

A 6-pin jack is located on the rear side of the device for connecting the PT 1000 series, contact thermometer or the contact plug. The electronics of the devices returns a test current that must flow via connector pins 3 and 5 for the heating plate to heat up.

Safety function:

If the test current is interrupted because of e.g. breakage of contact thermometer or falling out of the cable plug, the heating cuts off.

Settings:

For detailed instructions for settings and limit values, please refer to the operating instructions of the device you are connecting.

The desired medium temperature can be adjusted on the contact thermometer. The required surface temperature of the heating plate can be selected with the rotating / pressing knob or button.

Adjusting the temperature of device to the maximum adjustable temperature will result in the fastest possible heating time. However, the medium temperature may fluctuate to values above the set-point temperature on the contact thermometer. By adjusting the rotating / pressing knob or button to approximately twice the set-point value of contact thermometer (with a setpoint of + 60 °C, the temperature of device would be set to + 120 °C), you will reach a good compromise between a fast heating time and over-shooting the set point. If you adjust the temperature of the device to exactly the set-point temperature, the medium will not reach the set-point temperature because some loss of the heat will always occur between the heating plate and the medium.

The maximum heating plate temperature is limited to the adjusted safety temperature limit in the event of a control circuit malfunction. (See "Adjusting the safety temperature limit")

The rated temperature is setting on the VT 4.

If you are using an VT 4, we recommend using the VT 4 in On/ Off mode (two- point- control) at higher medium temperatures.

Error messages

Version VMS- D:

The LED displays (C and D) will start to flash and display a code when a fault is present. The fault codes and troubleshooting procedure are as follows:

Error code	LED	Cause	Solution
E2	G	Temperature inside device too high	- Switch off device and allow it cool down. - Check that the ventilation slots on the bottom are not blocked and that cooling air can circulate freely.
-5-	G	Safety circuit is triggered.	- Reduce the temperature setpoint or reset the safety circuit where necessary.
E2	D	No pulse to the reading fork (faulty reading fork, faulty motor or seized motor)	- Switch off the unit. - Trace the external cause blocking the stirring drive. - If you cannot find and visible cause, clean the unit and return it for repair.
E6	G	Not put contact plug Break and/or short-circuit of the contact thermometers	- Put contact plugs - Exchange the contact thermometer
Other codes			- Switch off the unit. - Clean and return for repair.

Version VMS- A:

The warning light (I) flashes to indicate that a fault is present. To remove a fault condition, proceed as follows:

- Switch off the unit and allow it to cool down.
- Trace and remove the cause of the fault. (Insert contact plug or contact thermometer)
- Switch on the unit.

If the fault persists, clean the unit and return it for repair.

Mounting support rod H 16 V

The H 16 V support rod can be screwed into the threading on the right behind the housing as follows:

- Remove the plastic cap and washer from the support rod and screw off the hex nut.
- Place the washer and hex nut on the beginning of the threading on the housing and hold them in place.
- Screw the support rod into the nut and threading of the housing and counter-tighten the hex nut.

Maintenance and cleaning

Version VMS- A or VMS- D:

The device is maintenance-free. It is only subject to the natural wear and tear of components and their statistical failure rate.

Cleaning

- For cleaning disconnect the mains plug!
- Use only cleaning agents which have been approved by VWR to clean VWR devices:
Water containing surfactant / isopropyl alcohol.
- Wear protective gloves during cleaning the devices.
- Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.
- Do not allow moisture to get into the device when cleaning.
- Before using another than the recommended method for cleaning or decontamination, the user must ascertain with VWR that this method does not destroy the device.

Spare parts order

When ordering spare parts, please give:

- device type.
- serial number, see type plate.
- position number and description of spare part, see www.vwr.com.

Repair

Please only send devices in for repair that have been cleaned and are free of materials which might present health hazards.

For repair, please request the "Safety Declaration (Decontamination Certificate)" from VWR or use the download printout of it from VWR website at www.vwr.com.

If your appliance requires repair, return it in its original packaging. Storage packaging is not sufficient when sending the device - also use appropriate transport packaging.

Version VMS- D

Information for care and maintenance of the heating plate with technical enamel coating:

The technical enamel makes the heating plate easier to care for and more resistant to acids and bases. Because of it, however, the heating plate is also more susceptible to extreme fluctuations in temperature and the force of impact. This can result in cracks forming or the coating flaking off.

Make certain that the bottom of the placing vessel is even, clean and dry. The bottom of the placing vessel must not have any sharp grooves, sides or edges. Remove residues of bases and immediately. We recommend most strongly that you clean the heating plate regularly.

Accessories

For accessories see www.vwr.com.

Technical Data

		VMS- A	VMS- D
Design voltage	VAC	220 - 240 ±10%	220 - 240 ±10%
or	VAC	115 ±10%	115 ±10%
or	VAC	100 ±10%	100 ±10%
Design frequency	Hz	50/60	50/60
Motor power consumption	W	15	15
Motor power output	W	2	2
Speed range	rpm	100 to 2000, infinitely variable setting	100 to 2000, infinitely variable setting
Adjustment of the speed control		rotary knob on the front	rotary knob on the front
Display of the speed control		scale 0 - 6	LED-display four digit (0 - 2000 rpm)
Motor		shaped pole motor	shaped pole motor
Max. stirring volume (water)	ltr	10	15
Heating plate		1.4301	1.4301 with enamel coating
Hating plate dimensions	mm	Ø 125	130 x 130
Heating power	W	400 (230 VAC)	500 (230 VAC)
	W	400 (115 VAC)	500 (115 VAC)
	W	360 (100 VAC)	450 (100 VAC)
Temperature range	°C	RT ... 320	RT ... 320
Temperature setting		rotary knob on the front	rotary knob on the front
Temperature display		scale 0 - 6	LED-display four digit (0°C ... +320°C)
Contact thermometer connection		-	DIN 12 878
Infinitely variable safe temperature	°C	100 ... 400	-
Fixed adjustment safe temperature	°C	400	-
Perm. ambient temperature	°C	+5 to +40	+5 to +40
Perm. humidity	%	80	80
Perm. on time	%	100	100
Screened for radio interference		under VDE 0871	under VDE 0871
Protection to DIN 60529		IP 21	IP 21
Contamination level		2	2
Overvoltage categorie		II	II
Protection class		1(protective earth)	1 (protective earth)
Fuse	A	10AT Id-Nr. 2755400	10AT Id-Nr. 2755400
Operation at a terrestrial altitude	m	max. 2000 above sea level	max. 2000 above sea level
Dimensions		168 x 220 x 105	168 x 220 x 105
Weight		2,4	2,4

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India

VWR Lab Products Private Limited
135/12, Brigade Towers, 2nd Floor
Front wing, Brigade Road,
Bengaluru, India – 560 025
Tel: +91-2522-647911/922 (Mumbai)
Tel: +91-80-41117125/26 (Bangalore)
Fax +91-80-41117120
E-mail: vwr_india@vwr.com

Ireland / Northern Ireland

VWR International Ltd / VWR
International (Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin, Dublin 15
Tel.: 01 88 22 222
Fax: 01 88 22 333
E-mail sales@ie.vwr.com

Italy

VWR International PBI S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: 02-3320311/02-487791
Fax: 02-332031307/02-40090010
E-mail: info@it.vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: 020 4808 400
Fax: 020 4808 480
E-mail: info@nl.vwr.com

New Zealand

VWR International LP
241 Bush Road
Albany
0632 New Zealand

Norway

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: 02290
Fax: 815 00 940
E-mail: info@no.vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: 058 32 38 210
Fax. 058 32 38 205
E-mail: labart@pl.vwr.com

Portugal

VWR International - Material de
Laboratório, Lda
Edifício Neopark
Av. Tomás Ribeiro, 43- 3 D
2790-221 Carnaxide
Tel.: 21 3600 770
Fax: 21 3600 798/9
E-mail: info@pt.vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel: +65 6505 0760
Fax: +65 6264 3780
E-mail: sales@sg.vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnologia 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: 902 222 897
Fax: 902 430 657
E-mail: info@es.vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: 08 621 34 00
Fax: 08 621 34 66
E-mail: info@se.vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: 044 745 13 13
Fax: 044 745 13 10
E-mail: info@ch.vwr.com

Turkey

Pro Lab Laboratuvar Teknolojileri Ltd.Şti.
a VWR International Company
Orta Mah. Cemal Gürsel Caddesi
Ördekcioglu İşmerkezi No.32/1
34896 Pendik - Istanbul
Tel.: +90216 598 2900
Fax: +90216 598 2907
Email: info@pro-lab.com.tr

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard
Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: 0800 22 33 44
Fax: 01455 55 85 86
E-mail: uksales@uk.vwr.com