BINDER

PURE AQUA SERVICE

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Water treatment system

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Mounting Instructions

Manuel de montage

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BINDER GmbH

Address

Tel.

Fax

Internet

E-mail

Service Hotline

Service Fax

Service E-mail

Service Hotline USA

Postfach 102,

D-78502 Tuttlingen

+49 7462 2005 0

+49 7462 2005 100

<http://www.binder-world.com>

info@binder-world.com

+49 7462 2005 555

+49 7462 2005 93 555

service@binder-world.com

+1 866 816 8191

Inhalt

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Content

1.	Safety	19
1.1.	Structure of the safety instructions.....	19
1.2.	General safety instructions on installing and operating BINDER Pure Aqua Service	20
1.3.	Material safety data sheet of the cartridge content.....	21
1.4.	Intended use	22
2.	Description of the equipment	22
3.	Scope of delivery, transportation, storage, and installation.....	22
3.1.	Scope of delivery.....	22
3.2.	Transport and storage.....	23
3.3.	Installation and ambient conditions.....	23
4.	Mounting.....	24
4.1.	Water connection of the cartridge	24
4.2.	Connection of the measuring equipment	25
4.3.	Connection to the climatic chamber.....	26
5.	Start up	28
6.	Indications on the measuring equipment	29
7.	Replacing the cartridge	30
8.	Service / contact	30
9.	Performance, lifetime	30
9.1.	Water hardness.....	30
9.2.	Water consumption	31
10.	Decommissioning.....	31
11.	Disposal.....	32
11.1.	Disposal of the transport packing.....	32
11.2.	Disposal of the cartridge	32
11.3.	Disposal of the measuring equipment and the mains unit in the Federal Republic of Germany.....	32
11.4.	Disposal of the measuring equipment and the mains unit in the member states of the EC except for the Federal Republic of Germany	33
11.5.	Disposal of the measuring equipment and the mains unit in non-member states of the EC.....	33

1. Safety

These mounting instructions are part of the scope of delivery. Always keep them at hand. To avoid injuries and damage, observe the safety instructions given with the mounting instructions.

	 CAUTION
	Failure to observe the safety instructions may result in serious injuries and unit damage. ➤ Observe the safety instructions in these mounting instructions ➤ Carefully read the complete mounting instructions

1.1. Structure of the safety instructions

In these mounting instructions, the following harmonized denominations and symbols indicate dangerous situations following the harmonization of ISO 3864-2 and ANSI Z535.6.

Signal word panel

Depending on the seriousness and probability of the consequences, dangers are identified with a signal word, the corresponding safety color, and if appropriate, the safety alert symbol.

 WARNING
Indicates an imminently hazardous situation that, if not avoided, will result in death or serious (irreversible) injury.

 CAUTION
Indicates a potentially hazardous situation which, if not avoided, may result in moderate or minor (reversible) injury.

CAUTION
Indicates a potentially hazardous situation which, if not avoided, may result in damage to the product and/or its functions or of a property in its proximity.

Safety alert symbol

Use of the safety alert symbol indicates **risk of injury**.

Observe all measures that are marked with the safety alert symbol in order to avoid death or injury.

Pictograms

Warning signs			
			
Electrical hazard	Danger of frost	Harmful substances	Pressurized device
			
Explosive substances			
Mandatory action signs		Prohibition signs	
			
Read instructions	Environment protection	Non-potable water	

Word message panel structure

Type / cause of hazard.

Possible consequences.

- Ø Instruction how to avoid the hazard: prohibition
- Instruction how to avoid the hazard: mandatory action

Observe all other notes and information not necessarily emphasized in the same way, in order to avoid disruptions which could result in direct or indirect injury or property damage.

Please observe the information given on the cartridge label.

1.2. General safety instructions on installing and operating BINDER Pure Aqua Service

NO liquid must pass by the cable connection into the inside of the measuring equipment.

	 WARNING
	Electrical hazard. Danger of death. Damage to the equipment. ➤ Make sure that the installation site is dry.

To avoid any possibility of water damage, ensure that there is a floor drain in the vicinity of where the instrument is located.

Water treated with BINDER Pure Aqua Service is not adapted as drinking water. Drinking the treated water can implicate loss of electrolytes.

	CAUTION
	Non-potable water. Health hazard. Ø Do NOT drink processed water.

The cartridge must not freeze since it was filled with water the first time.

	CAUTION
	Danger of frost. Damage to the cartridge. Ø Do NOT expose the cartridge to temperatures below +3°C / 37.4°F.

The connected cartridge is a pressurized device. Do not open. Secure against tipping over. The content may irritate eyes and skin.

 	CAUTION
	Opening by force or damage to the cartridge. Risk of injury. Irritation of skin and eyes. Damage to the cartridge and its environment. Ø Do NOT open. Ø Prevent from tipping over.

After using the equipment, close off the water supply at the freshwater supply.

1.3. Material safety data sheet of the cartridge content

	You can order the complete material safety data sheet of the cartridge content at BINDER service.
--	--

The following indications refer to the cartridge content. They are only relevant if the content emerges, p.ex due to damage of the cartridge.

Extract from the material safety data sheet of the cartridge content

- Content: Functionalized styrene / divinylbenzene copolymer: Polystyrene sulphonic acid and polyvinyl benzyl dimethyl amine or polystyrene sulphonic acid and polyvinyl benzyl trimethyl ammonium hydroxide.
- Avoid eye and skin contact. Keep away from food and beverages. Eye contact: rinse with plenty of water and seek medical attention. Contact with skin: wash with plenty of water. Ingestion: drink 200-300 ml (half pint) water and seek medical attention
- Do not discharge the cartridge content into drains.

1.4. Intended use

The water treatment system BINDER Pure Aqua Service serves for tap water treatment, preparing water suitable for operating BINDER climatic chambers. Using the water treatment system BINDER Pure Aqua Service is recommended with a local water hardness exceeding 142.8 mg/l (ppm) = 10 English (Clark) degrees = 8.32° grains per U.S. gallon = 1.4285 mmol/l.

Water treated with the water treatment system BINDER Pure Aqua Service is not suitable for analytical and / or medical purpose. It is no drinking water.

2. Description of the equipment

The water treatment system BINDER Pure Aqua Service consists of a reusable einer wiederverwendbaren measuring equipment to assess water quality, an exchangeable cartridge for water treatment, and the necessary connecting elements.

It serves for tap water treatment.

Dimensions of cartridge: height 67 cm / 26.4 in, width 22 cm / 8.7 in

Weight Pure Aqua Service with transport packing (state of delivery): 16 kg / 35.3 lb

Weight of cartridge not filled with water: 13 kg / 28.7 lb

Weight of cartridge filled with water: 19 kg / 41.9 lb

Power consumption of measuring equipment: 0.24 W up to 0.40 W

3. Scope of delivery, transportation, storage, and installation

3.1. Scope of delivery

The water treatment system BINDER Pure Aqua Service (Art. no. 8012-0625) consists of:

- Disposable cartridge (Art. no. 6011-0077)
- Measuring equipment mounted to T-piece, with mains unit 110V-240V and socket adapter
- Mounting instructions (Art. no. 7001-0159)
- Water hose for processed water 3m, screwing, and 2 gaskets
- Water drain for pressure compensation when replacing the cartridge

Figure 1: Scope of delivery BINDER Pure Aqua Service

3.2. Transport and storage

Permissible ambient temperature range during transport and storage:

- Cartridge not filled with water (state of delivery): -10°C / 14°F up to $+40^{\circ}\text{C}$ / 104°F .
- Cartridge filled with water: $+3^{\circ}\text{C}$ / 37.4°F up to $+40^{\circ}\text{C}$ / 104°F . Recommended storage temperature: max. 10°C / 50°F .

The cartridge must not freeze since it was filled with water the first time.

	CAUTION
	Danger of frost. Damage to the cartridge. Ø Do NOT expose the cartridge to temperatures below $+3^{\circ}\text{C}$ / 37.4°F.

Permissible ambient humidity: 70 % r.H. max., non-condensing

Intermediate storage in a closed and dry room.

3.3. Installation and ambient conditions

Set up the cartridge upright on an even surface. The connected cartridge is a pressurized device. Secure against tipping over. The content may irritate eyes and skin.

 	 CAUTION
	Opening by force or damage to the cartridge. Risk of injury. Irritation of skin and eyes. Damage to the cartridge and its environment. Ø Do NOT open. Ø Prevent from tipping over.

Permissible ambient temperature range during operation: $+3^{\circ}\text{C}$ / 37.4°F up to $+40^{\circ}\text{C}$ / 104°F (climatic chamber: $+18^{\circ}\text{C}$ / 64.4°F to $+32^{\circ}\text{C}$ / 89.6°F .)

- Permissible ambient humidity: 70 % r.H. max., non-condensing.
- Installation height: max. 2000 m above sea level.

To avoid any possibility of water damage, ensure that there is a floor drain in the vicinity of where the instrument is located!

To completely isolate the unit from the mains supply, you must disconnect the mains plug. Install the unit in a way that the mains unit is easily accessible and can be easily disconnected in case of danger.

The unit must not be installed and operated in potentially explosive areas.

	 WARNING
	Explosion hazard. Danger of death. Ø Do NOT operate the unit in potentially explosive areas. ➤ KEEP explosive dust or air-solvent mixtures AWAY from the vicinity of the unit.

4. Mounting

4.1. Water connection of the cartridge

Freshwater temperature: not below +5°C and not above 40°C.

Supply pressure: 1 bar up to 10 bar

Admissible substances: Max admissible chlorine content: 0.1 mg/l.

- Unscrew the yellow covers from the cartridge. You can use them for transport of a used the cartridge.
- Connect the water drain to the inlet of the cartridge. The inlet is marked with an arrow in the direction of the water flow.

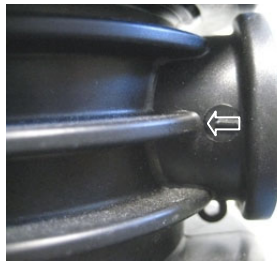
Figure 2: Cartridge inlet

Figure 3: Cartridge inlet with connected water drain

- Connect the freshwater hose delivered with the climatic chamber (R3/4" connecting piece, brass) to the water drain.

Figure 4: Cartridge with connected freshwater hose

- Connect the other end of the freshwater hose (R3/4" connecting piece, brass) to a lockable freshwater tap.

4.2. Connection of the measuring equipment

- Connect the measuring equipment which is mounted to a T-piece to the outlet of the cartridge and fix it with the dark grey R3/4" union nut. The outlet is marked with an arrow in the direction of the water flow.

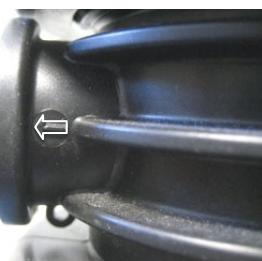
Figure 5: Cartridge outlet

- Connect the mains unit to the measuring equipment and plug it into a socket with a distribution voltage between 110V and 240V / 50/60 Hz.

Figure 6: Measuring equipment connected to cartridge and mains unit

No liquid must pass by the cable connection into the inside of the measuring equipment.

	WARNING Electrical hazard. Danger of death. Damage to the equipment. ➤ Make sure that the installation site is dry.
---	---

4.3. Connection to the climatic chamber

- Connect the hose for purified water (R3/4" plastic connecting piece) to the R3/4 in male thread of the measuring equipment. Insert one of the two gaskets delivered with the water hose. Conduct the hose to the climatic chamber.

Figure 7: Delivered gaskets at the hose for purified water (state of delivery)

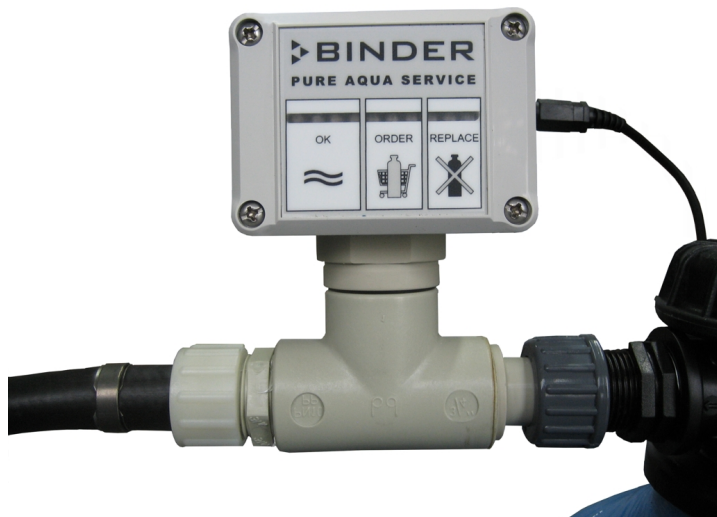
Figure 8: Measuring equipment connected to cartridge and mains unit with hose for purified water

- Connect the hose (R3/4" plastic connecting piece) to the freshwater connection at the rear of the climatic chamber. Insert one of the two gaskets delivered with the water hose. Observe the indications on water connection given in the chamber's operating manual.

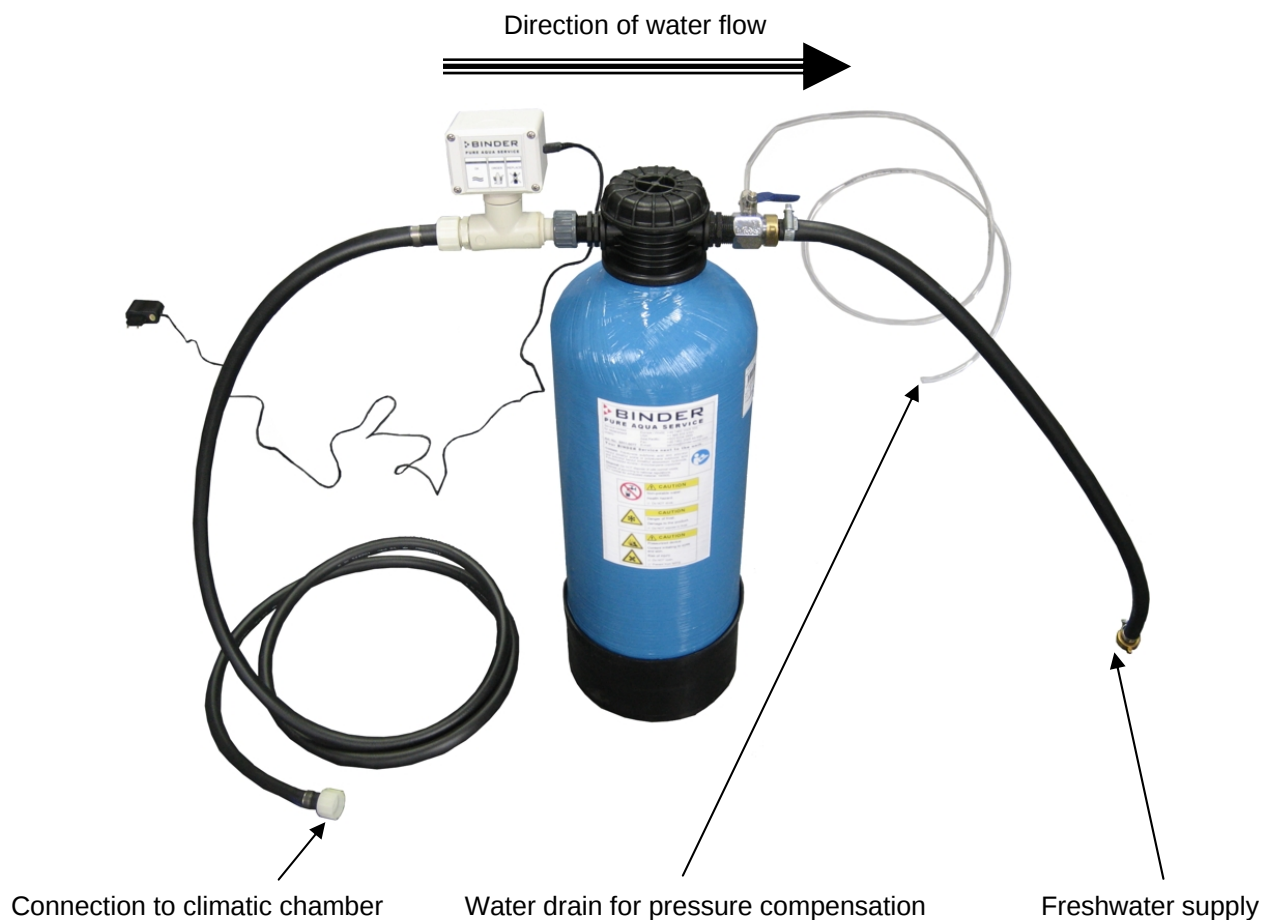
Figure 9: Disposable cartridge with measuring equipment and hoses, readily connected

Figure 10: KBF with connected water treatment system BINDER Pure Aqua Service

5. Start up

The cartridge must stand upright during use. Prevent it from tipping over.

- Close the water drain

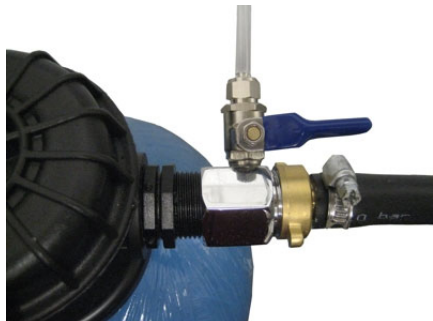
Figure 11: Closed water drain
(to operate the cartridge)

Figure 12: Open water drain
(to replace the cartridge)

- Slowly turn on the freshwater supply tap.

The water treatment system BINDER Pure Aqua Service is now ready for use.

With a new cartridge, the first green lamp on the measuring equipment lights up (chap. 6).

No liquid must pass by the cable connection into the inside of the measuring equipment.

	WARNING Electrical hazard. Danger of death. Damage to the equipment. ➤ Make sure that the installation site is dry.
--	---

Water treated with BINDER Pure Aqua Service is not adapted as drinking water. Drinking the treated water can implicate loss of electrolytes.

	CAUTION Non-potable water. Health hazard. Ø Do NOT drink processed water.
--	---

The connected cartridge is a pressurized device. Do not open. Secure against tipping over. The content may irritate eyes and skin.

 	CAUTION Opening by force or damage to the cartridge. Risk of injury. Irritation of skin and eyes. Damage to the cartridge and its environment. Ø Do NOT open. Ø Prevent from tipping over.
--	---

After using the equipment, close off the water supply to the measuring equipment at the freshwater tap.

6. Indications on the measuring equipment

	Lighting of the green lamps Very good water quality
	Additional lighting of the yellow lamps Very good water quality. Order a replacement cartridge Art. No. 6011-0077 at BINDER service (chap. 8).
	Additional lighting of the red lamps Still adequate water quality. Replace the cartridge. When all 3 red lamps are lit, the adequate water quality is no longer assured.

Note: With a longer duration of use, the time after which the next lamp lights up diminishes.

7. Replacing the cartridge

- Close the freshwater supply tap
- Introduce the free end of the tube, which is connected to the water drain into a receptacle of a volume of at least 0.5 liters / 0.018 cu.ft.
- Open the water drain. Due to the overpressure inside the cartridge, approx. 100 ml water will run off.
- Close the water drain. Remove all connections from the cartridge.
- Close the cartridge inlet and outlet for transport with the yellow covers. You can now dispose of the cartridge (chap. 11.2).

8. Service / contact

Order number for replacement cartridge: **6011-0077**

Contact BINDER Service for orders:

Europe / World: +49 7462 2005 555

USA: +1 866 816 8191

Asia Pacific: +603 6204 2855

Fax: +49 7462 2005 93 555

E-mail: service@binder-world.com

9. Performance, lifetime

The measuring equipment to assess the water quality is reusable.

The lifetime of the disposable cartridge depends on the water quality (water hardness) and the water consumption. We actually estimate the average lifetime at 1 to 1.5 years.

If the freshwater contains salts due to a preceding water treatment installation or additions of chlorine, the lifetime of the cartridge may be reduced. Max. admissible chlorine content: 0.1 mg/l. Desalinization plants for sea water have no influence.

9.1. Water hardness

Water hardness (freshwater)	Cartridge sufficient for processing the following quantity of water
178.5 mg/l (ppm) = 12.5 English (Clark) degrees = 10.4° grains per U.S. gallon = 1.785 mmol/l	approx. 2000 l
356 mg/l (ppm) = 25 English (Clark) degrees = 20.8° grains per U.S. gallon = 3.57 mmol/l	approx. 1000 l
712 mg/l (ppm) = 50 English (Clark) degrees = 41.6° grains per U.S. gallon = 7.14 mmol/l	approx. 500 l

International conversion table for water hardness:

Concentration of alkaline earths ions mmol/l	Concentration of calcium carbonate (CaCO ₃) mg/l = ppm *	Concentration of calcium oxide (CaO) mg/l	German degrees °dH	French degrees °f	English degrees / Clark degrees °E / °Clark	Concentration of CaCO ₃ grain per U.S. gallon gpg *
1	100,0	56.2	5.62	10.0	7.026	5.845
0,017	1.78	1	0.10	0.178	0.125	0.104
0,1785	17.85	10	1	1.785	1.25	1.04
0,10	10.0	5.6	0.56	1	0.700	0.582
0,14	14.2	8.0	0.8	1.428	1	0.832
0,17	17.1	9.6	0.96	1.71	1.20	1
1.428	142.8	80.0	8.0	14.28	10	8.32
1,785	178,5	100,0	10,0	17,85	12,5	10,4
3,57	356	200,0	20,0	35,6	25	20,8
5,355	534	300,0	30,0	53,4	37,5	31,2
7,14	712	400,0	40,0	71,2	50	41,6

* Both types figure as "American degrees".

9.2. Water consumption

Standard values of water consumption

KBF 240, KBF-ICH 240, KBF-LQC 240, KBWF 240:

- with set-points 25°C / 77°F and 60% r.H.: approx. 1 l per day (without opening the door)
- with set-points 40°C / 104°F and 75% r.H.: approx. 2 l per day (without opening the door)

MKF 240:

- with set-points 25°C / 77°F and 60% r.H.: approx. 1 l per day (without opening the door)
- with set-points 55°C / 131°F and 60% r.H.: approx. 2 l per day (without opening the door)

With humidity cycles, water consumption can increase up to a daily amount of 20 l.

10. Decommissioning

Close the water supply at the freshwater tap. Disconnect the measuring equipment from the mains. Remove the connections at the freshwater supply and at the climatic chamber.

- Temporal decommissioning: See indications for appropriate storage, chap. 3.2
- Final decommissioning: Dispose of the unit as described in chap. 11.

When taking into operation the system again, please observe the indications given in chap. 5.

11. Disposal

11.1. Disposal of the transport packing

Packing element	Material	Disposal
Transport box with metal clamps	Cardboard	Paper recycling
	Metal	Metal recycling
Foamed plastic stuffing	PE foam	Plastic recycling
Top cover	Cardboard	Paper recycling
Edge protection	Styropor®	Plastic recycling
Protection of doors and racks	PE foam	Plastic recycling
Bag for operating manual	PE foil	Plastic recycling
Insulating air cushion foil	PE foil	Plastic recycling

If recycling is impossible, all packing parts can also be disposed of with normal waste.

11.2. Disposal of the cartridge

- Do not dispose of the cartridge with normal waste.
- Dispose the content of the cartridge according to the local and national regulations on a disposal site or in compliance with the synthetic material methods. Dispose to an authorized waste collection point. A controlled combustion is possible. Do not discharge into drains or the environment. In state of delivery (cartridge not filled with water), the material is not classified as hazardous waste (exception: UK).
- EU number for exhausted or saturated material used for the preparation of drinking water or water for industrial use is 19 09 05
- EU number for exhausted or saturated material used in waste water treatment plants not otherwise specified is 19 08 06
- In UK, surplus product should be declared a 'Special Waste'. Refer to the 'Control of Pollution (Special Waste) Regulations 1980 - SI 1709'

11.3. Disposal of the measuring equipment and the mains unit in the Federal Republic of Germany

According to directive 2002/96/EC of the European Parliament and of the Council on waste electrical and electronic equipment (WEEE), BINDER devices are classified as "monitoring and control instruments" (category 9) only intended for professional use". They must not be disposed of at public collecting points.

The measuring equipment and the mains unit bear the symbol for the marking of electrical and electronic equipment manufactured / placed on the market in the EC after 13 August 2005 and be disposed of in separate collection according to the directive 2002/96/EC on waste electrical and electronic equipment (WEEE) and German national law for electrical and electronic equipment (Elektro- und Elektronikgerätegesetz, ElektroG). WEEE marking: crossed-out wheeled bin with solid bar under. A significant part of the materials must be recycled in order to protect the environment.


At the end of the device's service life, have the device disposed of according to the German national law for electrical and electronic equipment (Elektro- und Elektronikgerätegesetz, ElektroG) from 23 March 2005, BGBl. I p. 762 or contact BINDER service who will organize taking back and disposal of the unit according to the German national law for electrical and electronic equipment (Elektro- und Elektronikgerätegesetz, ElektroG) from 23 March 2005, BGBl. I p. 762.

	CAUTION
	Violation against existing law. ⊘ Do NOT dispose of BINDER devices at public collecting points. ➤ Have the device disposed of professionally at a recycling company which is certified according to the German national law for electrical and electronic equipment (Elektro- und Elektronikgerätegesetz, ElektroG) from 23 March 2005, BGBl. I p. 762. <i>or</i> ➤ Instruct BINDER service to dispose of the device. The general terms of payment and delivery of BINDER GmbH apply, which were valid at the time of purchasing the unit.

11.4. Disposal of the measuring equipment and the mains unit in the member states of the EC except for the Federal Republic of Germany

According to directive 2002/96/EC of the European Parliament and of the Council on waste electrical and electronic equipment (WEEE), BINDER devices are classified as "monitoring and control instruments" (category 9) only intended for professional use". They must not be disposed of at public collecting points.

The measuring equipment and the mains unit bear the symbol for the marking of electrical and electronic equipment manufactured / placed on the market in the EC after 13 August 2005 and be disposed of in separate collection according to the directive 2002/96/EC on waste electrical and electronic equipment (WEEE). WEEE marking: crossed-out wheeled bin with solid bar under.

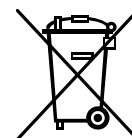At the end of the device's service life, notify the distributor who sold you the device, who will take back and dispose of the unit according to the directive 2002/96/EC of 27 January 2003 on waste electrical and electronic equipment (WEEE).

	CAUTION
	Violation against existing law. ⊘ Do NOT dispose of BINDER devices at public collecting points. ➤ Have the device disposed of professionally at a recycling company which is certified according to conversion of the directive 2002/96/EC into national law. <i>or</i> ➤ Instruct the distributor who sold you the device to dispose of it. The agreements apply that were reached with the distributor when purchasing the unit (e.g. his general terms of payment and delivery). ➤ If your distributor is not able to take back and dispose of the unit, please contact BINDER service.

11.5. Disposal of the measuring equipment and the mains unit in non-member states of the EC

	CAUTION
	Alteration of the environment. ➤ For final decommissioning and disposal of the climatic chamber, please contact BINDER service. ➤ Observe the statutory regulations for appropriate, environmentally friendly disposal.

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