# VWR<sup>®</sup> Thermometer

## INSTRUCTION MANUAL

### North American Catalog Numbers:

**12777-838, 89094-782, 89094-768, 89094-770**

### European Catalogue Numbers:

**620-1582, 620-1583, 620-1584, 620-1585**

Version: 1

Issued: 15<sup>th</sup> May 2014

# Legal Address of Manufacturer

## United States

VWR International  
Radnor Corporate Center  
Building One, Suite 200  
100 Matsonford Road, P.O. Box 6660  
Radnor, PA 19087  
800-932-5000  
<http://www.vwr.com>

## Europe

VWR International bvba  
Researchpark Haasrode 2020  
Geldenaaksebaan 464  
B-3001 Leuven  
+ 32 16 385011  
<http://be.vwr.com>

Country of origin – China

## Product Specifications

Range: -50 to 70°C (-58 to 158°F)

Resolution: 0.1°

Cat. No	Accuracy (°C)	Probe Type	Probe Length	Cable Length
VWRI620-1582	±0.5	Bottle	63.5 mm	3 m
VWRI620-1583	±0.5	Sensor	19 mm	3 m
VWRI620-1584	±1.0	Bottle	63.5 mm	3 m
VWRI620-1585	±1.0	Sensor	19 mm	3 m

## Operation - English

### TEMPERATURE DISPLAY MODES

**Ambient Temperature Display Mode (IN):** Indicated by “IN” appearing on the display, displays the current ambient temperature and either the ambient temperature memory (MIN/MAX) or the ambient temperature alarm limits (LO/HI).

### Probe Temperature Display Mode (OUT):

Indicated by “OUT” appearing on the display, displays the current probe temperature and either the probe temperature memory (MIN/MAX) or the probe temperature alarm limits (LO/HI). To change between ambient (IN) and probe (OUT) temperature display modes, press the IN/OUT button located on the back of the unit. To change between displaying the memory (MIN/MAX) and the alarm limits (HI/LO), press the MODE button.

### DISPLAYING °F OR °C

To display the temperature readings in Fahrenheit or Celsius, press the °C/°F button located on the back of the unit.

## MINIMUM AND MAXIMUM MEMORY

There are four points that are automatically recorded into memory:

- Minimum (MIN) Ambient Temperature Achieved
- Maximum (MAX) Ambient Temperature Achieved
- Minimum (MIN) Probe Temperature Achieved
- Maximum (MAX) Probe Temperature Achieved

All four memories are independent of each other.

**Minimum and maximum temperature memories are NOT programmable.**

The minimum temperature recorded into memory is the minimum temperature achieved since the last time the memory was cleared. The maximum temperature recorded into memory is the maximum temperature achieved since the last time the memory was cleared. The minimum and maximum temperature memories are maintained over the period since the memory was cleared.

Minimum and maximum memories are recorded for both ambient and probe temperatures regardless of the display mode.

## CLEARING THE PROBE (OUT) TEMPERATURE MEMORY

1. Select the Probe Temperature Display Mode by pressing the IN/OUT button located on the back of the unit until "OUT" appears on the display.
2. Press the MODE button until "MIN" and "MAX" appears on the display.
3. Press the MIN-RESET button to clear the minimum temperature memory.
4. Press the MAX-RESET button to clear the maximum temperature memory.

## CLEARING THE AMBIENT (IN) TEMPERATURE MEMORY

1. Select the Ambient Temperature Display Mode by pressing the IN/OUT button located on the back of the unit until "IN" appears on the display.
2. Press the MODE button until "MIN" and "MAX" appears on the display.
3. Press the MIN-RESET button to clear the minimum temperature memory.
4. Press the MAX-RESET button to clear the maximum temperature memory.

## ALARMS

Both ambient (IN) and probe (OUT) temperature alarm limits may be set in 1° increments.

With the alarm enabled (🔔), the unit will sound an alarm when the temperature being displayed is outside the alarm limits that have been set for that display mode (equal to or lower than the low alarm set point, or equal to or greater than the high alarm set point). Example: If the unit is displaying the ambient temperature (IN) and the probe temperature (OUT) is outside the probe temperature alarm limits, no alarm will sound.

## SETTING PROBE (OUT) TEMPERATURE ALARM LIMITS

1. Select the Probe Temperature Display Mode by pressing the IN/OUT button located on the back of the unit until "OUT" appears on the display.
2. Press MODE button until "LO" and "HI" appears.
3. To set the low temperature alarm limit (LO), press the LO ALM SET button to advance the display until the desired low temperature is displayed.

4. To set the high temperature alarm limit (HI), press the HI ALM SET button to advance the display until the desired high temperature is displayed.

### SETTING AMBIENT (IN) TEMPERATURE ALARM LIMITS

1. Select the Ambient Temperature Display Mode by pressing the IN/OUT button located on the back of the unit until "IN" appears on the display.
2. Press MODE button until "LO" and "HI" appears.
3. To set the low temperature alarm limit (LO), press the LO ALM SET button to advance the display until the desired low temperature is displayed.
4. To set the high temperature alarm limit (HI), press the HI ALM SET button to advance the display until the desired high temperature is displayed.

### ENABLE/DISABLE ALARMS

To enable the alarm to sound when the temperature being displayed is outside the alarm limits, press the  button located on the back of the unit until "" appears on the display. Each successive press of the  button will enable () or disable the alarm.

### ALARM SOUNDING

With the alarm enabled () , the unit will sound an alarm when the temperature being displayed is outside the alarm limits that have been set for that display mode (equal to or lower than the low alarm set point, or equal to or greater than the high alarm set point). If the alarm is sounding based on the low alarm limit, "LO" will flash on the display. If the alarm is sounding based on the high alarm limit, "HI" will flash on the display. The unit will continue to alarm and flash the corresponding "LO" or "HI" until the temperature being displayed is within the alarm limits. Once the temperature being displayed is within the alarm limits, the alarm will stop sounding and the corresponding "LO" or "HI" will no longer flash.

To silence an alarm manually, press any button. When manually silenced, the corresponding "LO" or "HI" will continue to flash on the display until the temperature is within the alarm limits. When enabled () , the alarm will only sound if the temperature being displayed is outside the alarm limits that have been set for that display mode (IN or OUT). Example: If the unit is displaying ambient temperature (IN) and the probe temperature (OUT) is outside the probe temperature alarm limits, **no alarm will sound**.

### RESET BUTTON

The RESET button located on the back of the unit may be used to simultaneously reset the memory and alarm limits for both the ambient (IN) and probe (OUT) sensor.

To reset the unit, use a pointed object to press the RESET button.

When the reset button is pressed, the following will occur:

- The display will show all segments for approximately 5 seconds.
- The ambient (IN) and probe (OUT) minimum/maximum memory will be cleared.
- The ambient (IN) and probe (OUT) alarm limits will be returned to the default value of 10°C (LO) and 30°C (HI).
- The unit will display the ambient (IN) temperature in °C.

## ON/OFF BUTTON

The ON/OFF button located on the back of the unit may be used to turn the unit off when the thermometer is not being used. The minimum/maximum memory and alarm limits are not cleared when the unit is turned off. To turn the unit off, press the ON/OFF button. Each successive press of the ON/OFF button will turn the unit on or off.

**No alarms will sound and no minimum/maximum memory will be recorded when the unit is off.**

## TROUBLESHOOTING

If this unit does not function properly for any reason, replace the battery with a new high-quality battery (see the “Battery Replacement” section). Low battery power can occasionally cause any number of “apparent” operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

## TRACEABLE® CERTIFICATE

To assure accuracy, the unit is supplied with an individually serial-numbered Traceable® Certificate provided from an ISO/IEC 17025 calibration laboratory accredited by A2LA. The recommended recalibration date is noted on the certificate that is included with the product. Annual recalibrations should occur thereafter.

## BATTERY REPLACEMENT

Erratic readings, a faint display, no display, or a battery symbol appearing on the display are all indications that the battery must be replaced. Open the battery compartment by depressing the tab located at the top of the battery cover and lifting it out. The battery cover is hinged at the bottom so that the battery compartment may be opened without completely removing the battery cover. Remove the exhausted battery and replace with a AAA alkaline battery. Make certain to insert the new battery with the proper polarity as indicated in the illustration in the battery compartment. Close the battery cover.

Replacing the battery has the same effect as pressing the RESET button. (See the “RESET BUTTON” section.)

## Technical service/Web Resources

Visit the VWR’s website at [www.vwr.com](http://www.vwr.com) for:

- Complete technical service contact information
- Access to VWR’s Online Catalogue, and information about accessories and related products
- Additional product information and special offers

**Contact us** For information or technical assistance contact your local VWR representative or visit.

[www.vwr.com](http://www.vwr.com).

## Warranty

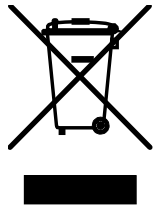
**VWR International** warrants that this product will be free from defects in material and workmanship for a period of two (2) years from manufacturing date. If a defect is present, VWR will, at its option, repair, replace, or refund the purchase price of this product at no charge to you, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear.

For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products. IT IS EXPRESSLY AGREED THAT THIS

WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

### **Equipment disposal**

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste. Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment. By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

### **Compliance with local laws and regulations**

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

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## Local VWR offices in Europe and Asia Pacific

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### India

VWR Lab Products Private Limited  
135/12, Brigade Towers, 2nd Floor  
Front wing, Brigade Road,  
Bengaluru, India - 560 025  
Tel: +91-2522-647911/922 (Mumbai)  
Tel: +91-80-41117125/26 (Bangalore)  
Fax +91-80-41117120  
E-mail: vwr\_india@vwr.com

### Ireland / Northern Ireland

VWR International Ltd / VWR International  
(Northern Ireland) Ltd  
Orion Business Campus  
Northwest Business Park  
Ballycoolin  
Dublin 15  
Tel.: 01 88 22 222  
Fax: 01 88 22 333  
E-mail sales@ie.vwr.com

### Italy

VWR International PBI S.r.l.  
Via San Giusto 85  
20153 Milano (MI)  
Tel.: 02-3320311/02-487791  
Fax: 02-332031307/02-40090010  
E-mail: info@it.vwr.com

### The Netherlands

VWR International B.V.  
Postbus 8198  
1005 AD Amsterdam  
Tel.: 020 4808 400  
Fax: 020 4808 480  
E-mail: info@nl.vwr.com

### New Zealand

VWR International LP  
241 Bush Road  
Albany  
0632 New Zealand

### Norway

VWR International AS  
Haavard Martinsens vei 30  
0978 Oslo  
Tel.: 02290  
Fax: 815 00 940  
E-mail: info@no.vwr.com

### Poland

VWR International Sp. z o.o.  
Limbowa 5  
80-175 Gdansk  
Tel.: 058 32 38 210  
Fax. 058 32 38 205  
E-mail: labart@pl.vwr.com

### Portugal

VWR International - Material de  
Laboratório, Lda  
Edifício Neopark  
Av. Tomás Ribeiro, 43- 3 D  
2790-221 Camaxide  
Tel.: 21 3600 770  
Fax: 21 3600 798/9  
E-mail: info@pt.vwr.com

### Singapore

VWR Singapore Pte Ltd  
18 Gul Drive  
Singapore 629468  
Tel: +65 6505 0760  
Fax: +65 6264 3780  
E-mail: sales@sg.vwr.com

### Spain

VWR International Eurolab S.L.  
C/ Tecnología 5-17  
A-7 Llinars Park  
08450 - Llinars del Vallès  
Barcelona  
Tel.: 902 222 897  
Fax: 902 430 657  
E-mail: info@es.vwr.com

### Sweden

VWR International AB  
Fagerstagatan 18a  
163 94 Stockholm  
Tel.: 08 621 34 00  
Fax: 08 621 34 66  
E-mail: info@se.vwr.com

### Switzerland

VWR International GmbH  
Lerzenstrasse 16/18  
8953 Dietikon  
Tel.: 044 745 13 13  
Fax: 044 745 13 10  
E-mail: info@ch.vwr.com

### Turkey

Pro Lab Laboratuvar Teknolojileri Ltd.Şti.  
a VWR International Company  
Orta Mah. Cemal Gürsel Caddesi  
Ördekioğlu İşmerkezi No.32/1  
34896 Pendik - Istanbul  
Tel.: +90216 598 2900  
Fax: +90216 598 2907  
Email: info@pro-lab.com.tr

### UK

VWR International Ltd  
Customer Service Centre  
Hunter Boulevard  
Magna Park Lutterworth  
Leicestershire  
LE17 4XN  
Tel.: 0800 22 33 44  
Fax: 01455 55 85 86  
E-mail: uksales@uk.vwr.com